

Setting up ImageScope on Mac

This application note discusses how to install ImageScope on a Mac.

Prerequisites

- **Intel Mac** – iMac, Mac Mini, Mac Book, Mac Book Pro, Mac Pro Tower
- Mac OS X “Tiger” 10.4.6 or higher running on Intel powered Mac
- Minimum memory requirement of 512 MB RAM, recommended 1 GIG Ram
- 30 MB available hard disk space for Parallels Desktop installation, plus sufficient space allocated for Windows XP operating system. 15 GIG of hard disk space recommended.
- **Parallels Desktop License** (see below).
- **Windows XP OS License and installation disk.**

Installing Parallels Desktop


- Download Parallels Desktop from www.parallels.com and obtain an activation key. You can download a 30 day evaluation copy prior to purchase, this is recommended as it provides you with the Parallels Tools.
- The instructions in this Application Note refer to build 3186 of Parallels desktop.
- Follow the instructions for installation in the Parallels Desktop for Mac Quick Start Guide found in the Parallels –Desktop folder on the desktop
- Follow the instructions in the Parallels Desktop for Mac Getting Started Guide up to Chapter 4 – Activating Parallels Desktop.


Parallels Desktop for Mac

Getting Started Guide




Installing a Virtual Machine


- Launch Parallels Desktop by double-clicking on the orange Parallels icon 
- Follow the instructions in the OS Installation Assistant.

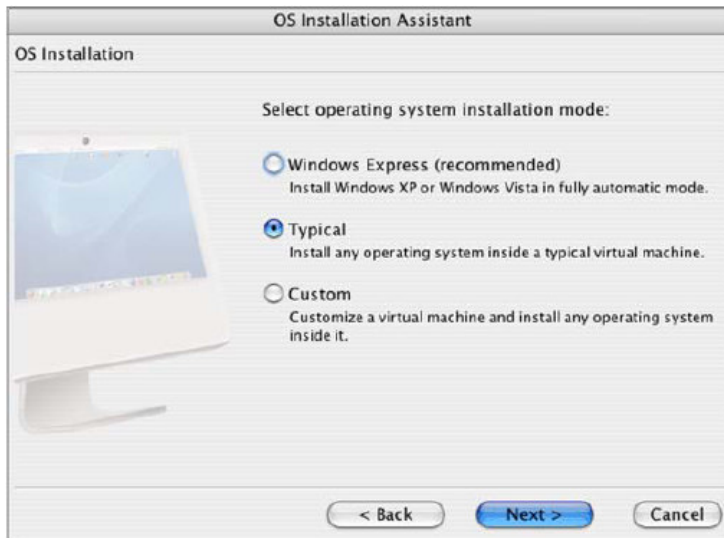

- Select Typical Windows OS Installation
- Choose Windows XP as the operating system. You will need an XP installation disk and product key (you will be asked for this later).

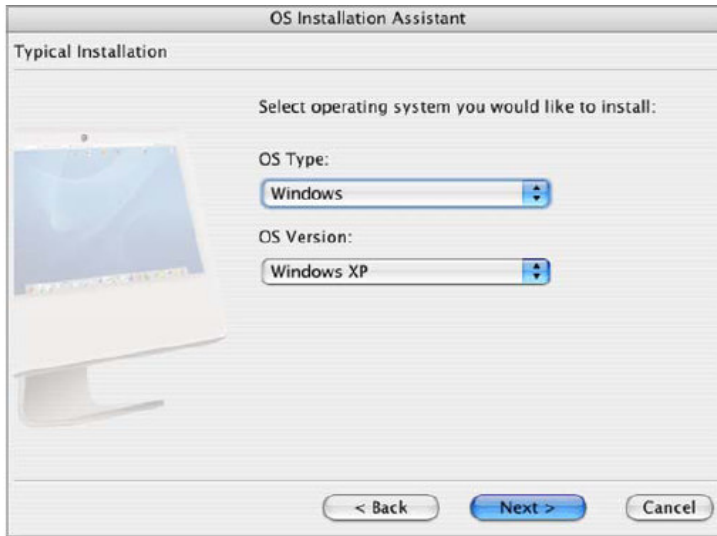

- Set the amount of RAM for the new virtual machine, but do not allocate any more than half of the physical RAM installed on your Mac.

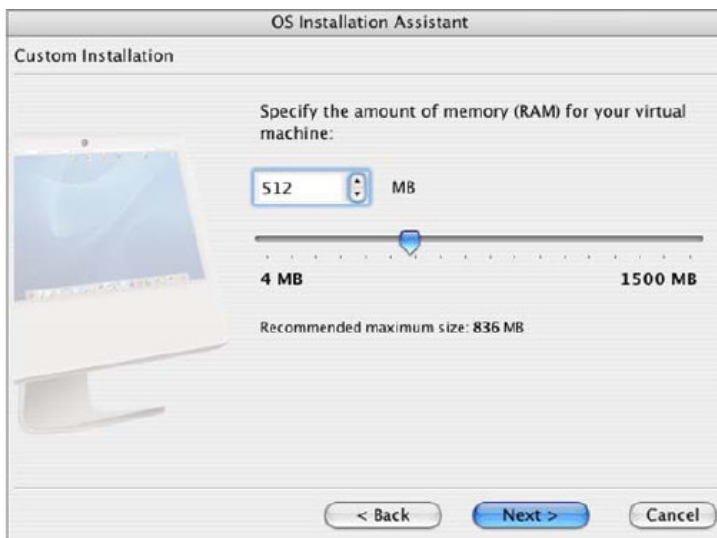

- Next choose whether you want to create a new virtual hard disk or not (refer to the Parallels for Mac User Guide, Chapter 4, for more details)



- Choose a name for your Virtual Machine.



- Put your XP disk in the CD/DVD Drive and follow the instructions.




- The XP installation will launch inside a new window and this will take some time to install (approximately 1 hour)
- Follow the on screen instructions and reference the instructions from Chapter 4 in the Parallels Desktop for Mac Getting Started Guide.
- When the installation has completed, install the Parallels Tools following the instructions from Chapter 4 (page64) in the Parallels Desktop for Mac Getting Started Guide.

Installing ImageScope

- Download ImageScope from www.aperio.com if you do not already have the ImageScope.exe file.
- Copy the ImageScope.exe installer onto a CD or flash drive.
- Open the XP Virtual Machine you created.
- Install ImageScope by clicking on the executable and following the instructions.

Opening Images in ImageScope

- Open the Virtual Machine you created.
- Images can be opened in several ways:
 - Open a browser in the Virtual Machine for example www.images2.scanscope.com and you will be able to open them by clicking on “Open with ImageScope”.
 - Create a shared folder of Images on the hard disk of your Mac using the Parallels Shared Folders tool, see Chapter 13 of the Parallels Desktop for Mac Getting Started Guide.

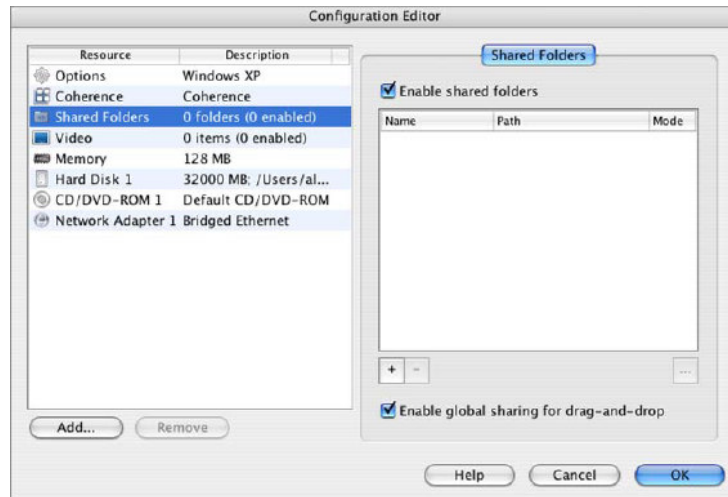

- Browse to this share inside the Virtual Machine in My Network Places and open Images by clicking on them.

IMPORTANT NOTE: ImageScope only works inside the Parallels Virtual Machine you created.

©Copyright 2007 Aperio Technologies, Inc.
Part Number/Revision: MAN-0019, Revision A
Date: 5th March, 2007

This document applies to software versions Release 8 and later.

All rights reserved. This document may not be copied in whole or in part or reproduced in any other media without the express written permission of Aperio Technologies, Inc. Please note that under copyright law, copying includes translation into another language.

User Resources

For the latest information on Aperio Technologies products and services, please visit the Aperio Technologies website at: <http://www.aperio.com>.

Disclaimers

This manual is not a substitute for the detailed operator training provided by Aperio Technologies, Inc., or for other advanced instruction. Aperio Technologies Field Representatives should be contacted immediately for assistance in the event of any instrument malfunction. Installation of hardware should only be performed by a certified Aperio Technologies Service Engineer.

ImageServer is intended for use with the svx file format (the native format for digital slides created by scanning glass slides with the ScanScope scanner). Educators will use Aperio software to view and modify digital slides in Composite WebSlide (CWS) format.

CAUTION

For research and education uses only, not for use in diagnostic procedures. This product has not been approved or cleared as a medical device by the U.S. Food and Drug Administration. The data and images obtained or viewed using this product are not intended for clinical or diagnostic use.

Trademarks and Patents

ScanScope is a registered trademark and ImageServer, TMA Lab, ImageScope, and Spectrum are trademarks of Aperio Technologies, Inc. All other trade names and trademarks are the property of their respective holders.

Aperio products are protected by U.S. Patents: 6,711,283; 6,917,696; 7,035,478; and 7,116,440; and licensed under one or more of the following U.S. Patents: 6,101,265; 6,272,235; 6,522,774; 6,775,402; 6,396,941; 6,674,881; 6,226,392; 6,404,906; 6,674,884; and 6,466,690.

Contact Information

Headquarters: Aperio Technologies, Inc.
1430 Vantage Court, Suite 106
Vista, CA 92081
United States

European Office: Aperio
3 The Sanctuary
Eden Office Park
Ham Green
Bristol BS20 0DD, UK

United States of America

Tel: +1 866-478-4111 (toll free)

Fax: +1 760-539-1116

Customer Service

Tel: +1 866-478-4111 (toll free)

Technical Support

Tel: +1 866-478-4111 (toll free)

Email: support@aperio.com

Europe

Tel: +44 (0) 1275 375123

Fax: +44(0) 1275 373501

Customer Service

Tel: +44 (0) 1275 375123

Technical Support

Tel: +44 (0) 1275 375123

Email: europesupport@aperio.com